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- 5) OECD (2008) Guidelines for the Testing of Chemicals Test No. 425: Acute Oral Toxicity: Up-and-Down Procedure. Available at:
<https://ntp.niehs.nih.gov/iccvm/suppdocs/feddocs/oeed/oeedtg425.pdf>
- 6) Halle, W. (2003) The registry of cytotoxicity: Toxicity testing in cell cultures to predict acute toxicity (LD₅₀) and to reduce testing in animals, *Alternatives to Laboratory Animals*, 31, 89-198.
- 7) NICEATM, NIEHS (2006) *In vitro* Cytotoxicity Test Methods for Estimating Acute Oral Systemic Toxicity BACKGROUND REVIEW DOCUMENT. Available at:
<https://ntp.niehs.nih.gov/pubhealth/evalatm/test-method-evaluations/acute-systemic-tox/in-vitro-validation/index.html>
- 8) ICCVAM, NICEATM (2006) ICCVAM TEST METHOD EVALUATION REPORT *In Vitro* Cytotoxicity Test Methods for Estimating Starting Doses for Acute Oral Systemic Toxicity Testing. Available at:https://ntp.niehs.nih.gov/iccvm/docs/acutetox_docs/brd_tmer/at-tmer-complete.pdf
- 9) OECD (2010) Guidance Document on using Cytotoxicity Tests to Estimate Starting Doses for Acute Oral Systemic Toxicity Tests. Series on Testing and Assessment No. 129. Available at:
<https://ntp.niehs.nih.gov/pubhealth/evalatm/test-method-evaluations/acute-systemic-tox/in-vitro-validation/index.html>
- 10) JaCVAM 急性毒性試験代替法評価委員会, 急性毒性試験代替法の第三者評価報告書 評価対象試験: *In vitro* 細胞毒性試験による急性経口毒性試験の初回投与量設定試験 (2011 年 1 月 14 日) Available at: https://www.jacvam.jp/files/list/07/07_01_C1.pdf
- 11) JaCVAM 評価会議, 単回投与毒性試験代替法の評価会議報告書 (2011 年 4 月 20 日). Available at: https://www.jacvam.jp/files/list/07/07_01_B1.pdf
- 12) ECVAM (2011) Follow-up study on the predictive capacity of the 3T3 Neutral Red Uptake cytotoxicity assay to correctly identify substances not classified for acute oral toxicity under the EU CLP system (LD₅₀>2000mg/kg) Final study report. Available at:
<https://www.jacvam.jp/list.html>
- 13) Prieto P., Cole, T., Curren R., b, Gibson R. M., Liebsch M., Raabe H., Tuomainen, A. M., Whelan, M. and Kinsner-Ovaskainen, A. (2013) Assessment of the predictive capacity of the 3T3 Neutral Red Uptake cytotoxicity test method to identify substances not classified for acute oral toxicity (LD₅₀>2000 mg/kg): results of an ECVAM validation study. *Regulatory Toxicology and Pharmacology*, 65(3) 344-365.
- 14) EURL ECVAM (2013) EURL ECVAM Recommendation on the 3T3 Neutral Red Uptake Cytotoxicity Assay for Acute Oral Toxicity. Available at:
<https://ec.europa.eu/jrc/en/publication/eur-scientific-and-technical-research-reports/eurl-ecvam-recommendation-3t3-neutral-red-uptake-cytotoxicity-assay-acute-oral-toxicity>
- 15) JaCVAM 急性毒性試験資料編纂委員会, 評価報告書 急性経口毒性を予測するための *In vitro* 細胞毒性試験 (2018 年 12 月 11 日). Available at:

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Concern approach in food safety assessment. Available at:

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- 1) National Institute for Occupational Safety and Health (1987) Registry of Toxic Effects of Chemical Substances 1985-86 Edition, Vol. 3, p.2535.
- 2) European Chemicals Agency, Information on chemicals, EC No. 202-494-5.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/1158/7/3/2>
- 3) Registry of Toxic Effects of Chemical Substances (RTECS) No. VJ2275000.
- 4) U.S. National Library of Medicine, ChemIDplus.
Available at: <https://chem.nlm.nih.gov/chemidplus/rn/50-99-7>
- 5) U.S. National Library of Medicine, ChemIDplus.
Available at: <https://chem.nlm.nih.gov/chemidplus/rn/57-50-1>
- 6) Cosmetic Ingredient Review (2015) Safety assessment of polysaccharide gums as used in cosmetics. Available at: <https://www.cir-safety.org/sites/default/files/plpogu092015final.pdf>
- 7) European Chemicals Agency, Information on chemicals, EC No. 201-788-0.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/13631/7/3/2>
- 8) European Chemicals Agency, Information on chemicals, EC No. 200-273-8.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/14143/7/3/2>
- 9) European Chemicals Agency, Information on chemicals, EC No. 200-811-1.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/13725/7/3/2>
- 10) European Chemicals Agency, Information on chemicals, EC No. 200-735-9.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/21618/7/3/2>
- 11) European Chemicals Agency, Information on chemicals, EC No. 200-291-6.
Available at: <https://echa.europa.eu/ro/registration-dossier/-/registered-dossier/14876/7/3/2>
- 12) Registry of Toxic Effects of Chemical Substances (RTECS) No. HA1600000.
- 13) European Chemicals Agency, Information on chemicals, EC No. 200-158-2.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/12475/7/3/2>
- 14) Registry of Toxic Effects of Chemical Substances (RTECS) No. MA2275100.
- 15) European Chemicals Agency, Information on chemicals, EC No. 200-293-7.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/12560/7/3/2>
- 16) Registry of Toxic Effects of Chemical Substances (RTECS) No. MB7600000.
- 17) European Chemicals Agency, Information on chemicals, EC No. 200-745-3.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/21605/7/3/2>
- 18) European Chemicals Agency, Information on chemicals, EC No. 200-798-2.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/20720/7/3/2>
- 19) European Chemicals Agency, Information on chemicals, EC No. 200-522-0.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/11897/7/3/2>
- 20) Registry of Toxic Effects of Chemical Substances (RTECS) No. OL5540000.

- 21) European Chemicals Agency, Information on chemicals, EC No. 211-519-9.
Available at: <https://echa.europa.eu/ro/registration-dossier/-/registered-dossier/12220/7/3/2>
- 22) European Chemicals Agency, Information on chemicals, EC No. 200-562-9.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/19745/7/3/2>
- 23) European Chemicals Agency, Information on chemicals, EC No. 200-568-1.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/16991/7/3/2>
- 24) European Chemicals Agency, Information on chemicals, EC No. 205-702-2.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/13104/7/3/2>
- 25) European Chemicals Agency, Information on chemicals, EC No. 200-274-3.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/5558/7/3/2>
- 26) European Chemicals Agency, Information on chemicals, EC No. 200-774-1.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/16986/7/3/2>
- 27) European Chemicals Agency, Information on chemicals, EC No. 200-795-6.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/10835/7/3/2>
- 28) U.S. National Library of Medicine, Pubchem.
Available at: <https://pubchem.ncbi.nlm.nih.gov/compound/Tyrosine#section=Acute-Effects>
- 29) European Chemicals Agency, Information on chemicals, EC No. 200-773-6.
Available at: <https://echa.europa.eu/registration-dossier/-/registered-dossier/14506/7/3/2>
- 30) 食品安全委員会（2004）食品健康影響評価の結果について，府食第 637 号，平成 16 年 6 月 10 日 . Available at: <https://www.fsc.go.jp/fscis/evaluationDocument/show/kya20040119160>
- 31) U.S. National Library of Medicine, ChemIDplus. Available at:
<https://chem.nlm.nih.gov/chemidplus/rn/107-21-1>
- 32) U.S. National Library of Medicine, ChemIDplus. Available at:
<https://chem.nlm.nih.gov/chemidplus/rn/107-88-0>
- 33) Cosmetic Ingredient Review (2018) Safety Assessment of alkane diols as used in cosmetics.
Available at: https://www.cir-safety.org/sites/default/files/alkane%20diols_0.pdf
- 34) European Chemicals Agency, Information on chemicals, EC No. 211-074-0. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/15109/7/3/235>
- 35) Cosmetic Ingredient Review (2015) Safety Assessment of glycerin as used in cosmetics.
Available at: https://www.cir-safety.org/sites/default/files/glycer_122014_FR.pdf
- 36) Cosmetic Ingredient Review (2010) Final report of the cosmetic ingredient review expert panel, Amended safety assessment of triethylene glycol and polyethylene glycols (PEGs) -4, -6, -7, -8, -9, -10, -12, -14, -16, -18, -20, -32, -33, -40, -45, -55, -60, -75, -80, -90, -100, -135, -150, -180, -200, -220, -240, -350, -400, -450, -500, -800, -2M, -5M, -7M, -9M, -14M, -20M, -23M, -25M, -45M, -65M, -90M, -115M, -160M and -180M and any PEGs ≥ 4 as used in cosmetics.
- 37) Cosmetic Ingredient Review (2012) Final report of the cosmetic ingredient review expert panel

- on the safety assessment of methyl acetate, International Journal of Toxicology, 31 (Supplement 1) 112S-136S.
- 38) European Chemicals Agency, Information on chemicals, EC No. 200-580-7. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/15549/7/3/2>
- 39) Registry of Toxic Effects of Chemical Substances (RTECS) No. ES5425000.
- 40) European Chemicals Agency, Information on chemicals, EC No. 203-532-3. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/13507/7/3/2>
- 41) U.S. National Library of Medicine, ChemIDplus. Available at:
<https://chem.nlm.nih.gov/chemidplus/rn/107-92-6>
- 42) 三菱化学テクノロジーリサーチ(2015)平成 27 年度報告毒物劇物指定のための有害性情報の収集・評価 物質名：ヘキサン酸. Available at:
<http://www.nihs.go.jp/law/dokugeki/hyouka/2015/142-62-1.pdf>
- 43) European Chemicals Agency, Information on chemicals, EC No. 205-550-7. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/14271/7/3/2>
- 44) Registry of Toxic Effects of Chemical Substances (RTECS) No. RH0175000.
- 45) European Chemicals Agency, Information on chemicals, EC No. 204-677-5. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/15370/7/3/2>
- 46) U.S. National Library of Medicine, ChemIDplus. Available at:
<https://chem.nlm.nih.gov/chemidplus/rn/124-07-2>
- 47) European Chemicals Agency, Information on chemicals, EC No. 203-931-2. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/13098/7/3/2>
- 48) European Chemicals Agency, Information on chemicals, EC No. 206-376-4. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/18512/7/3/2>
- 49) European Chemicals Agency, Information on chemicals, EC No. 205-582-1. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/15262/7/3/2>
- 50) European Chemicals Agency, Information on chemicals, EC No. 200-312-9. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/15218/7/3/2>
- 51) Registry of Toxic Effects of Chemical Substances (RTECS) No. PC1400000.
- 52) Registry of Toxic Effects of Chemical Substances (RTECS) No. KQ6300000.
- 53) Registry of Toxic Effects of Chemical Substances (RTECS) No. NT8050000.
- 54) Registry of Toxic Effects of Chemical Substances (RTECS) No. EO1400000.
- 55) European Chemicals Agency, Information on chemicals, EC No. 200-751-6. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/15322/7/3/2>
- 56) Registry of Toxic Effects of Chemical Substances (RTECS) No. EL5425000.
- 57) European Chemicals Agency, Information on chemicals, EC No. 204-633-5. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/13936/7/3/2/>

- 58) Registry of Toxic Effects of Chemical Substances (RTECS) No. MQ4025000.
- 59) U.S. National Library of Medicine, ChemIDplus. Available at:
<https://chem.nlm.nih.gov/chemidplus/rn/111-27-3>
- 60) European Chemicals Agency, Information on chemicals, EC No. 203-852-3. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/13265/7/3/2>
- 61) U.S. National Library of Medicine, ChemIDplus. Available at:
<https://chem.nlm.nih.gov/chemidplus/rn/111-70-6>
- 62) European Chemicals Agency, Information on chemicals, EC No. 203-897-9. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/5921/7/3/2>
- 63) European Chemicals Agency, Information on chemicals, EC No. 203-917-6. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/15210/7/3/2>
- 64) European Chemicals Agency, Information on chemicals, EC No. 203-982-0. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/15424/7/3/2>
- 65) European Chemicals Agency, Information on chemicals, EC No. 253-149-0. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/15942/7/3/2>
- 66) European Chemicals Agency, Information on chemicals, EC No. 203-751-4. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/16077/7/3/2>
- 67) Cosmetic Ingredient Review (2000) Final report, Safety assessment of glyceryl laurate, glyceryl laurate SE, glyceryl laurate/oleate, glyceryl adipate, glyceryl alginate, glyceryl arachidate, glyceryl arachidonate, glyceryl behenate, glyceryl caprate, glyceryl caprylate, glyceryl caprylate/caprate, glyceryl citrate/lactate/linoleate/oleate, glyceryl cocoate, glyceryl collagenate, glyceryl erucate, glyceryl hydrogenated rosinatate, glyceryl hydrogenated soyate, glyceryl hydroxystearate, glyceryl isopalmitate, glyceryl isostearate, glyceryl isostearate/myristate, glyceryl isostearates, glyceryl lanolate, glyceryl linoleate, glyceryl linolenate, glyceryl montanate, glyceryl myristate, glyceryl isotridecanoate/stearate/adipate, glyceryl oleate SE, glyceryl oleate/elaidate, glyceryl palmitate, glyceryl palmitate/stearate, glyceryl palmitoleate, glyceryl pentadecanoate, glyceryl polyacrylate, glyceryl rosinatate, glyceryl sesquioleate, glyceryl/sorbitol oleate/hydroxystearate, glyceryl stearate/acetate, glyceryl stearate/maleate, glyceryl tallowate, glyceryl thiopropionate, and glyceryl undecylenate.
- 68) Registry of Toxic Effects of Chemical Substances (RTECS) No. AH5425000.
- 69) Cosmetic Ingredient Review (1984) Final report on the safety assessment of candelilla wax, carnauba wax, Japan wax, and beeswax, Journal of the American College of Toxicology, 3 (3) 1-41.
- 70) Cosmetic Ingredient Review (2008) Safety Assessment of Simmondsia Chinensis (Jojoba) Seed Oil, Simmondsia Chinensis (Jojoba) Seed Wax, Hydrogenated Jojoba Oil, Hydrolyzed Jojoba Esters, Isomerized Jojoba Oil, Jojoba Esters, Simmondsia Chinensis (Jojoba) Butter, Jojoba

Alcohol, and Synthetic Jojoba Oil.

- 71) Cosmetic Ingredient Review (1980) Final report of the safety assessment for acetylated lanolin alcohol and related compounds, *Journal of Environmental Pathology and Toxicology* 4(4), 63-92.
- 72) European Chemicals Agency, Information on chemicals, EC No. 210-647-2. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/13795/7/3/2>
- 73) Cosmetic Ingredient Review (2018) Amended safety assessment of triglycerides as used in cosmetics.
- 74) U.S. National Library of Medicine, ChemIDplus. Available at:
<https://chem.nlm.nih.gov/chemidplus/rn/109-66-0>
- 75) Registry of Toxic Effects of Chemical Substances (RTECS) No. RZ9450000.
- 76) Registry of Toxic Effects of Chemical Substances (RTECS) No. MN9275000.
- 77) Cosmetic Ingredient Review (2012) Safety assessment of isoparaffins as used in cosmetics, *International Journal of Toxicology*, 31(Supplement 3) 269S-295S.
- 78) Japan Existing Chemical Database, Single dose toxicity test of 2,2,4,4,6,8,8-heptamethylnonane in rats. Available at: https://dra4.nihs.go.jp/mhlw_data/home/paper/paper4390-04-9a.html
- 79) Cosmetic Ingredient Review (1982) Final report on the safety assessment of squalane and squalene, *Journal of the American College of Toxicology*, 1 (2) 36-56.
- 80) Cosmetic Ingredient Review (2020) Safety assessment of polyene group as used in cosmetics, *International Journal of Toxicology*, 39 (Supplement 2) 59S-90S.
- 81) U.S. National Library of Medicine, ChemIDplus. Available at:
<https://chem.nlm.nih.gov/chemidplus/rn/63148-62-9>
- 82) European Chemicals Agency, Information on chemicals, EC No. 208-762-8. Available at:
<https://echa.europa.eu/registration-dossier/-/registered-dossier/15811/7/3/2>